



KERN & Sohn GmbH

Ziegelei 1

D-72336 Balingen

E-Mail: info@kern-sohn.com

Phone: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Installation Instructions

Foot switch

KERN KXS-A03

Version 1.0

08/2014

GB

KXS-A03-IA-e-1410



KERN KXS-A03

Version 1.0 08/2014

Installation Instructions

Foot switch





For KERN KXS-TM.


1 Scope of delivery

- Foot switch

2 Assembly

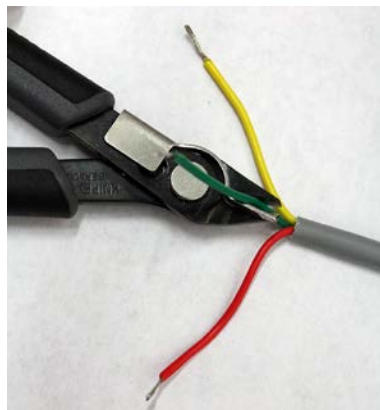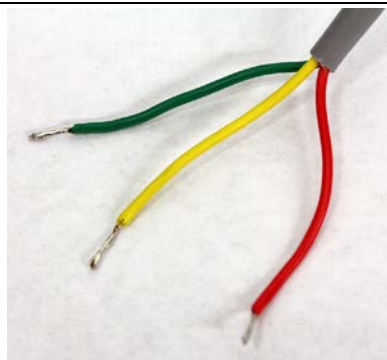



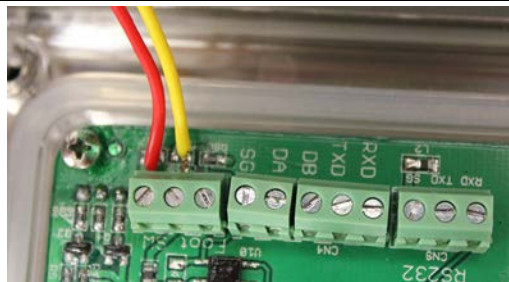




Cut off the green cable not necessary using pliers.




Connect the red and the yellow cable according to illustration.

3 Operation

The foot switch is used to send the weighing values to a printer or a PC. Alternatively can also be tared and the balance be reset to zero.

- ⇒ Start balance by pressing 
- ⇒ In the weighing mode press  and  at the same time, "01 FnC" is displayed
- ⇒ Press , „FnC 00“ is displayed.
- ⇒ Select „FnC 11“ using  and .
- ⇒ Acknowledge by .

FnC 11	<ul style="list-style-type: none">• ZERO• TARE• PRINT
---------------	---

- ⇒ Select the required function and acknowledge with 
- ⇒ In the weighing mode press foot switch, the selected function will be carried out